



FIBRE EXTRACTION SYSTEM F-6P Cellulose Analyser



Food Analysis

RELIABLE, VERSATILE AND ECONOMIC
FIBRE EXTRACTION SYSTEM FOR
EFFICIENT CRUDE AND DETERGENT
FIBRE ANALYSES

Fibre extraction system

Our fibre extraction system is specially optimized for the determination of Crude, Acid and Neutral fibre together with other fractions such as cellulose, hemicellulose

and lignin according to the universally known **Weende** and **Van Soest** methods.

Its flexibility to analyze a wide range of samples with subsequent or individual extraction, including boiling and filtration steps, makes our hot fibre extractor suitable for multiple settings and applications.


F-6P is designed to guarantee accurate and precise results in compliance with the international standardized methods such as **AOAC**, **AACC** and **ISO**

MAIN FIELDS OF APPLICATION

FOOD AND FEED INDUSTRY



Forage

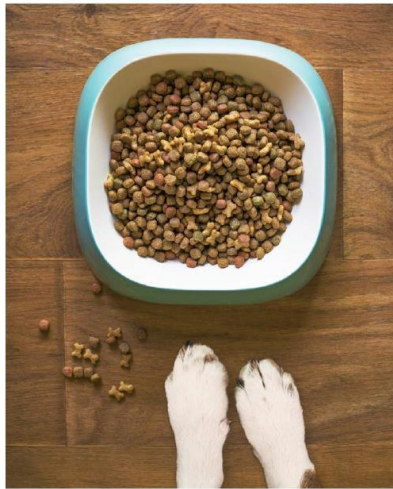

Pet food



Cereals

Office / Ofis

Email : export@bastak.com.tr
website: www.bastak.com.tr

Tel : +90 312 395 6 787
Fax: +90 312 395 6 788

İvedik OSB 1354.Cad 1387.Sk No:19
Yenimahalle - Ankara TURKEY

Factory / Fabrika

Dağyaka mahallesi 2010 .Cd.No:23
Kazan Ankara TURKEY

Tel : +90 312 811 2 811
Fax: +90 312 811 2 812

Email : info@bastak.com.tr
Website: www.bastak.com.tr



FEATURES

HIGHLY EFFICIENT

Simultaneous single or sequential hot extraction of 6 six samples under identical conditions including boiling and filtration.

TEMPERATURE CONTROL

Control of heat generation by a manual regulator. Guarantees rapid temperature rise until reagent boiling and temperature maintenance at lower levels as required.

EXEMPT FROM SAMPLE LOSS

Specialized heat resistant Pyrex® crucibles with filters are used for the extraction and filtration of components, preventing any sample loss as the crucibles can be used as sample vessels during extraction, weighing, drying and incineration.

REAGENT DISCHARGE

During the extraction the reagent can be discharged to the drain or collected for posterior fibre fraction analyses.

QUALITY AND SAFETY GUARANTEE

Easy to clean and corrosion resistant external frame made of stainless steel grade AISI-304 with epoxy coating. All the electrical and mechanical elements are fully protected.



Determination of multiple components.



Integrated peristaltic pump for a faster residue extraction.



Extraction and filtration without sample transfers.



High reproducibility of conditions and results.



According to standardized analysis procedures.



Personalized application support.



Integrated air pump for breaking compact clumps during filtration.



Easy to use.



Powerful quartz heater controlled by a regulator that homogeneously heats samples.

PRESSURE CONTRIBUTION

F-6P has an integrated peristaltic pump for negative pressure to discharge reagents and an air pump for positive pressure to break compact clumps during filtration.

Additionally, F-6P has an available positive pressure connection to EF-6P to break compact clumps during cold extractions, guaranteeing optimal working conditions while using both equipment.

VERSATILE USE

Multiple fibre portions are measurable across a wide variety of sample types. In each extraction phase, samples can be dried and quantified.

EFFICIENT EXTRACTION

Included peristaltic pump that allows the separation and elimination of reagents and residues quickly and efficiently, guaranteeing reliable and contamination-free analyses.


ACCORDING TO STANDARDIZED METHODS

Fibre content analysis with F-6P extractors are performed in accordance to official methods described by international entities such as AOAC, AACC and ISO in order to guarantee accurate results.



Fibre Extraction System

GENERAL OVERVIEW OF VAN SOEST METHOD FOR FIBRE COMPONENTS ISOLATION

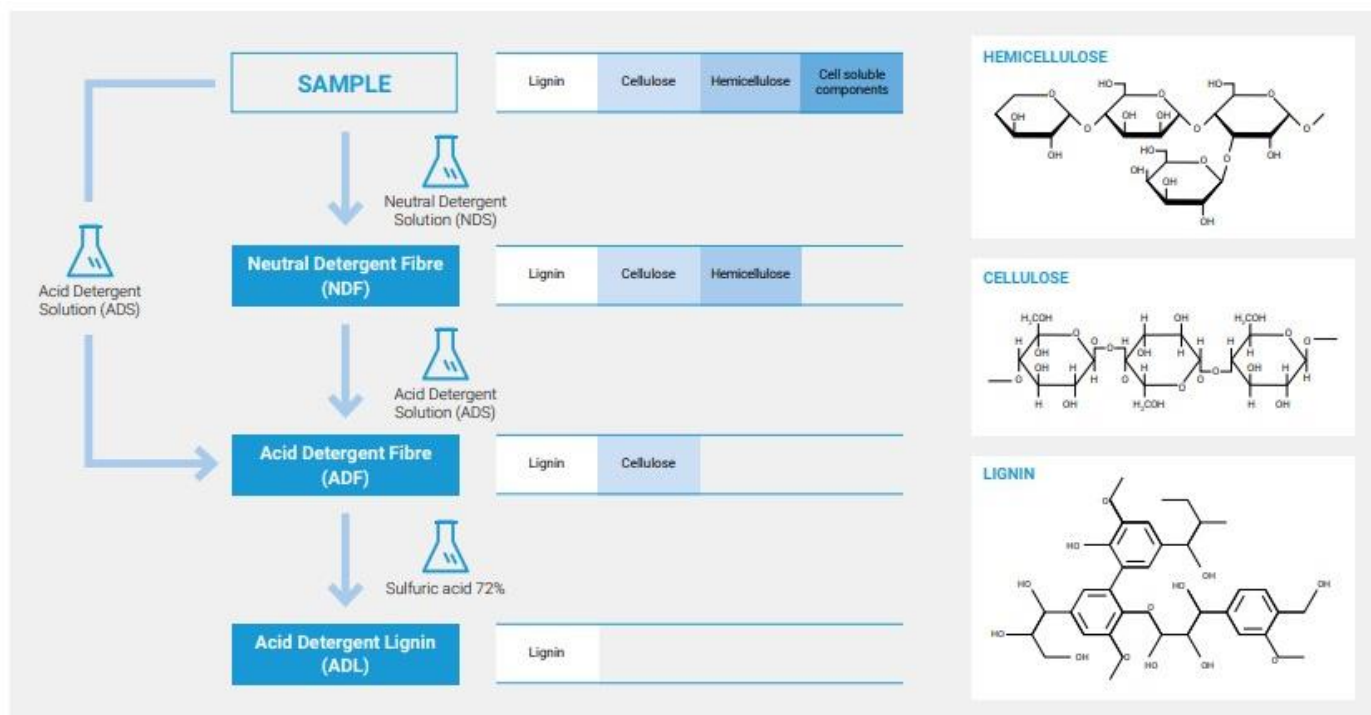

GENERAL OVERVIEW OF WEENDE METHOD FOR CRUDE FIBRE DETERMINATION



CRUDE FIBRE

Traditional crude fibre extraction, also known as the **Weende** method, is commonly used to estimate the quality of foods of plant origin on the premise that it constitutes their least digestible fraction, especially in monogastrics forage.

The analysis consists of a subsequent extraction with hot acid (1.25% H₂SO₄) and alkaline (1.25% KOH) solutions which removes protein, some hemi-cellulose and lignin.

Applications examples:
EN ISO 6865
AOAC 978.10

DETERGENT FIBRE

Detergent fibre is widely used to estimate the energy intake for ruminants. The analysis is done following the **Van Soest** method which is based on the principle that fibre can be further divided into less digestible fractions:

Acid Detergent Fibre (ADF)

Cell wall portions made up of cellulose and lignin. Indicates the amount of forage an animal can digest.

Applications examples:
EN ISO 13906
AOAC 973.18

Neutral Detergent Fibre (NDF)

Total cell wall including ADF and hemicelluloses. Indicates the amount of forage an animal can consume.

Applications examples:
ISO 16472
AOAC 2002:04

Acid Detergent Lignin (ADL)

Lignin fraction of ADF.

Applications examples:
EN ISO 13906
AOAC 973.18

Office/ Ofis

Email : export@bastak.com.tr
website: www.bastak.com.tr

Tel : +90 312 395 6 787
Fax: +90 312 395 6 788

İvedik OSB 1354.Cad 1387.Sk No:19
Yenimahalle Ankara TURKEY

Factory / Fabrika

Dağyaka mahallesi 2010 .Cd.No:23
Kazan Ankara TURKEY

Tel : +90 312 811 2 811
Fax: +90 312 811 2 812

Email : info@bastak.com.tr
Website: www.bastak.com.tr



MAIN FEATURES OF SOLUTION FOR FIBRE EXTRACTIONS

DURABLE EXTERNAL FRAME

- AISI-304 stainless steel
- Epoxy coating
- Easy to clean
- Corrosion resistant

WATER REFRIGERATION CIRCUIT

INTEGRATED AIR PUMP TO BREAK COMPACT CLUMPS (POSITIVE PRESSURE)

ADJUSTABLE TEMPERATURE REGULATION

PRESSURE CONNECTION TO F-6P TO BREAK COMPACT CLUMPS (POSITIVE PRESSURE)

6 SIMULTANEOUS EXTRACTION PLACES

DURABLE EXTERNAL FRAME

- AISI-304 stainless steel
- Epoxy coating
- Easy to clean
- Corrosion resistant

6 SIMULTANEOUS EXTRACTION PLACES

HEATING LID FOR A MORE HOMOGENEOUS HEAT DISTRIBUTION

INDIVIDUAL 3 POSITION VALVES

- Drain
- Closed
- Pressure support

INDIVIDUAL 3 POSITION VALVES

- Drain
- Closed
- Pressure support

INTEGRATED PERISTALTIC PUMP TO RECOVER SOLVENT (NEGATIVE PRESSURE)

INTEGRATED PERISTALTIC PUMP FOR RESIDUE EXTRACTION (NEGATIVE PRESSURE)

Office/ Ofis

Email : export@bastak.com.tr
website: www.bastak.com.tr

Tel : +90 312 395 6 787
Fax: +90 312 395 6 788

İvedik OSB 1354.Cad 1387.Sk No:19
Yenimahalle Ankara TURKEY


Factory / Fabrika

Dağyaka mahallesi 2010 .Cd.No:23
Kazan Ankara TURKEY

Tel : +90 312 811 2 811
Fax: +90 312 811 2 812

Email : info@bastak.com.tr
Website: www.bastak.com.tr

EXTRACTION STEPS FOR F-6P

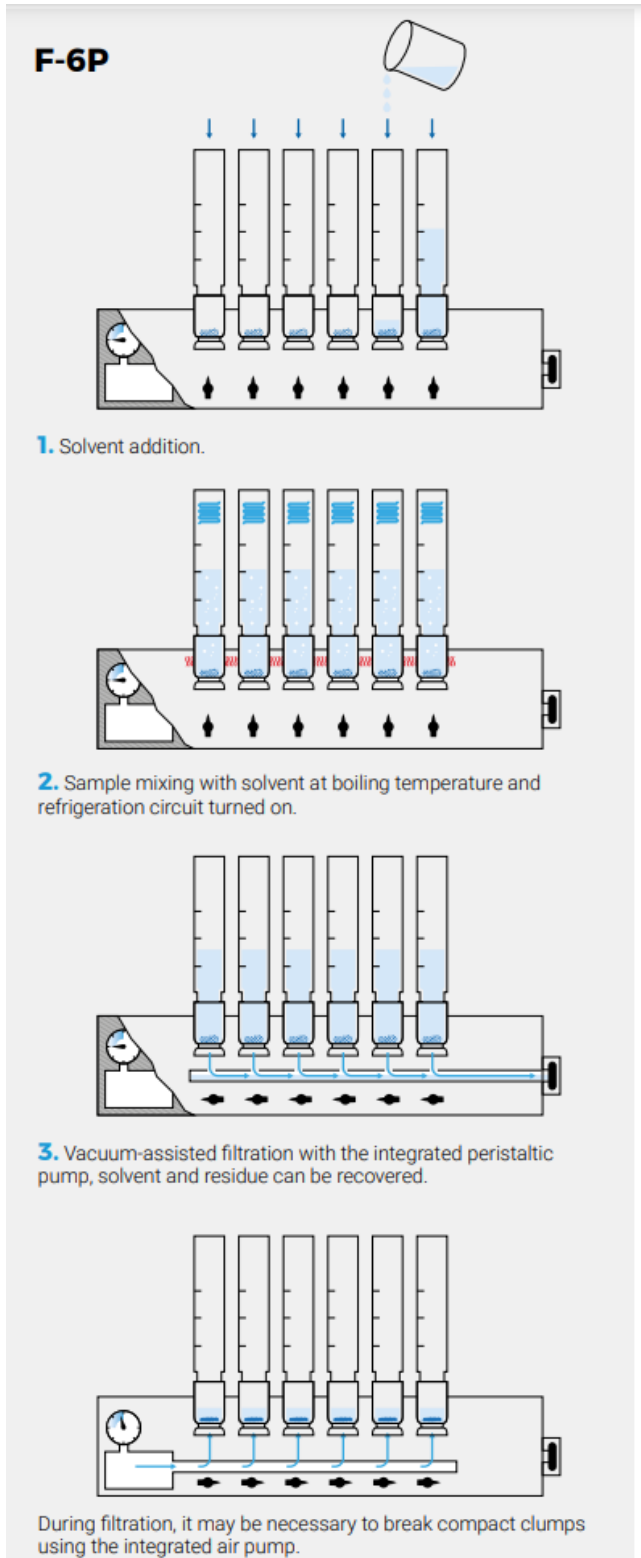

RECOMMENDED APPLICATIONS

- Determination of crude fibre content according to Weende.
- Determination of Acid Detergent Fibre (ADF) and Neutral Detergent Fibre (NDF) according to Van Soest.
- Determination of Acid Detergent Lignin (ADL).

TECHNICAL DESCRIPTION

- Simultaneous single or sequential hot or cold extraction of 6 six samples under identical conditions including boiling, rinsing and filtration.
- Integrated system without sample transfer or sample loss, samples can be dried or weighed on every stage via heat resistant crucibles.
- Heating by a quartz heater controlled by a manual regulator.
- Built-in water refrigeration circuit for hot extractions.
- Built-in air pump for positive pressure to break compact clumps during filtration, activated by an independent switch.
- Integrated peristaltic pump for negative pressure to drain solvent faster or optional fibre fraction collection, activated by an independent switch.
- Pyrex® crucibles with a nominal porosity of 40-90µm.
- Control of extraction steps by 3-position valves (Closed, Drain and Positive pressure).
- Gaskets and connection hoses made of Viton® compatible with various reagents including acid and base solutions.
- Easy-to-clean and corrosion resistant external frame made of AISI-304 stainless steel painted with epoxy resin.
- All electric and mechanical elements are conveniently protected.

SUPPLIED WITH THE FOLLOWING COMPONENTS:

- Crucibles set of 6 units.
- Crucibles rack.
- Tong for an individual manipulation of crucibles.
- Tong for a simultaneous manipulation of 6 crucibles.
- Heater lid.
- Several connection hoses.
- Several connection hoses clamps.



TECHNICAL SUMMARY OF F-6 P

	General info	Extractor general classification	Semiautomatic
	Extraction places	6	
	Standards compliance	AOAC, AACC, ISO	
	Dimensions L x D x H mm	724 x 330 x 580	
	Weight Kg	41	
	Power W	1250	
	Voltage V	230 V (115 V optional)	
	Frequency Hz	50/60	
	Recommended applications	Determination of Crude fibre content according to Weende	✓
	Determination of Acid Detergent Fibre (ADF) and Neutral Detergent Fibre (NDF) according to Van Soest	✓	
	Determination of Acid Detergent Lignin (ADL)	✓	
	Materials	Crucibles	Pyrex® glass
	Crucibles gaskets	Viton®	
	Boiling vessels	Borosilicate 3.3	
	Tubing	Silicone and Viton®	
	External housing	AISI-304 stainless steel painted with epoxy resin	
	Technological features	Quartz heater	✓
	Air pump (positive pressure to break compact clumps)	✓	
	Peristaltic pump (negative pressure for reagent discharge)	✓	
	Control panel	Temperature regulation	Switch + Rotary regulator
	Activation of positive pressure support to break compact clumps	Switch + Individual valves	
	Activation of negative pressure support to drain or recover reagents	Switch + Individual valves	
	Performance for fibre analysis	Sample capacity/batch units	6
	Sample capacity/day units	36	
	Sample quantity per crucible g	0,5 - 3	
	Precision (relative) %	± 0,1	
	Measuring range %	0,1 - 100	
	Reproducibility %	1 - 30	
	Functions gained with the accessory cold fat extraction system EF-6P	Preliminary degreasing before hot fibre extraction with F-6P	✓
	Cold fat extraction with acetone or other non-polar organic solvents	✓	



Complete batch handling items included

6 GLASS CRUCIBLES OF P2 POROSITY	✓
1 CRUCIBLES RACK	✓
1 TONG FOR AN INDIVIDUAL MANIPULATION OF CRUCIBLES	✓
1 TONG FOR A SIMULTANEOUS MANIPULATION OF 6 CRUCIBLES	✓
1 HEATER LID	✓
SEVERAL CONNECTION HOSES	✓
SEVERAL CONNECTION HOSES CLAMPS	✓



Accessories

CRUCIBLES SET

Reference	CR-P2
Dimensions Ø x H mm	34 x 60
Material	Pyrex® glass
Porosity grade	P2
Nominal porosity µm	40-90
Quantity units	6



COLD FAT EXTRACTION SYSTEM

Reference	EF-6P
External dimensions L x D x H mm	715 x 320 x 285
Power W	30
Voltage* V	230
Weight Kg	15
Frequency Hz	50/60
Analysis time min	40
Number of sample positions	6
Compatible crucibles dimensions Ø x H mm	34 x 60





Dimensions and performance



Reference	F-6P	EF-6P
External dimensions L x D x H mm	724 x 330 x 580	724 x 320 x 285
Power W	1280	30
Voltage* V	230	230
Weight Kg	41	15
Frequence Hz	50/60	50/60
Sample capacity	Batch units	6
	Day units	36
Glass crucible dimensions Ø x H mm	34 x 60	34 x 60

*Also available with a voltage of 115 V.

International standardized methods

F-6P hot fibre extractors are fabricated guaranteeing compliance with a variety of international standards such as AOAC, AACC and ISO.

Main fields of application

FOOD AND FEED INDUSTRY



FORAGE



PET FOOD



CEREALS

Regulations

Our F-6P fibre extraction system is designed to comply with the strictest international directives and standards, including the following:

- **EN-61010-1** Safety requirements for electrical equipment for measurement, control and laboratory use. **Part 1: General requirements.**
- **EN-61010-2-010 Part 2-010** Particular requirements for laboratory equipment for the heating of materials.
- **EN-61326** Electrical equipment for measurement, control and laboratory use. EMC Requirements.
- **2014/35/UE** Low voltage.
- **2014/30/UE** Electromagnetic compatibility.

When you work with BASTAK you receive;

- ✓ Free technical consultancy for tailor made solutions,
- ✓ High quality instruments with free webinars and on-line trainings,
- ✓ Fast delivery of instruments, spare parts and consumables,
- ✓ Remote technical support (Audio and Video) six (6) days a week,
- ✓ Economical spare parts and consumables available in 10 years,
- ✓ Free factory visits, live tests and demos,
- ✓ On-Site Warranty and On-Site After Sales Service (for >20.000 USD orders)

Office/ Ofis

Email : export@bastak.com.tr
website: www.bastak.com.tr

Tel : +90 312 395 6 787
Fax: +90 312 395 6 788

İvedik OSB 1354.Cad 1387.Sk No:19
Yenimahalle Ankara TURKEY

Factory / Fabrika

Dağyaka mahallesi 2010 .Cd.No:23
Kazan Ankara TURKEY

Tel : +90 312 811 2 811
Fax: +90 312 811 2 812

Email : info@bastak.com.tr
Website: www.bastak.com.tr



“More Than 23 Years of Experience ...”

Bastak Laboratory

57 Types of Device, 35 Types of Additive



Youtube > Bastak Expert

<https://www.youtube.com/watch?v=ktEz2PHdDQg&t=10s>

Bastak Factory - 7000 m2

20.000 instruments in 142 countries



Youtube > Bastak Factory

<https://www.youtube.com/watch?v=UbtWWhEbLy4>



You can find all Catalogues and Competitor Comparison Tables in this link:

https://drive.google.com/drive/folders/1_4aXqljmlSG0Gi7c0KeMuPAGWqvi65mT



Sensitive
Nutrition



High
Efficiency



Monoblock
Study Body



Standard
Grinding



Personal
Security

**The information in this document is subject to change without any prior notice or obligation on the part of the manufacturer and should not be construed as a commitment by the manufacturer. The manufacturer assumes no responsibility for any errors that may appear in this document.*

Office / Ofis

Email : export@bastak.com.tr
website: www.bastak.com.tr

Tel : +90 312 395 6 787
Fax: +90 312 395 6 788

İvedik OSB 1354.Cad 1387.Sk No:19
Yenimahalle Ankara TURKEY

Factory / Fabrika

Dağyaka mahallesi 2010 .Cd.No:23
Kazan Ankara TURKEY

Tel : +90 312 811 2 811
Fax: +90 312 811 2 812

Email : info@bastak.com.tr
Website: www.bastak.com.tr